



Precipio, Inc.

High Resolution Melting Protocol

High Resolution Melting of ICE COLD-PCR Products using the
ABI7500 Fast System

Table of Contents

| | |
|--|---|
| Principles of High Resolution Melting | 2 |
| High Resolution Melting of the ICE COLD-PCR products | 2 |
| Amplification Setup | 3 |
| ABI7500 Fast Instrument Setup | 4 |
| Review and Analysis of the HRM Melt Curve | 5 |
| Mutation Confirmation of HRM Positive Samples | 7 |
| Technical Support/Questions | 7 |

Copyright © 2018 by Precipio, Inc.

All rights reserved. This Instruction manual or any portion thereof may not be reproduced in any form, or transmitted outside of the recipient's organization in any form by any means — electronic, mechanical, photocopy, recording, or otherwise — or used in any manner whatsoever without the express written permission of the Company.