



Precipio, Inc.

Extraction Protocol

Extraction of DNA from plasma using the NextPrep-Mag™
cfDNA Isolation Kit

Table of Contents

Notes for this Assay	2
Reagents and Equipment Required but not Supplied	2
Workplace Preparation Suggestions	3
Plasma DNA Extraction Protocol	3

Copyright © 2018 by Precipio, Inc.

All rights reserved. This Instruction manual or any portion thereof may not be reproduced in any form, or transmitted outside of the recipient's organization in any form by any means — electronic, mechanical, photocopy, recording, or otherwise — or used in any manner whatsoever without the express written permission of the Company